



Compact, Inexpensive and Made for Small Scale Work

Whether you work in a research or a clinical laboratory, the compact **CAPPWash Microplate Washer** is the ideal washing tool for small scale immuno assays. Fitting into even the tightest budget, **CAPP Wash** is a perfect alternative to dependence on department washers for washing microplates and strips.



Features:

- CAPP Wash Microplate Washer performs uniformly and efficiently every time
- Washing is simple. Simply lower the tubes into the well, then push the knob to fill and release the knob to empty.
 Repeat for the desired number of washings and move on to the next row
- Matches all well types. The washer can be set to the desired depth very easily by simply adjusting the two finger screws found on either side of the tubes
- The CAPP Wash is available in 8-channel, 12-channel and 16-channel versions, as well as a 6-channel Tube Wash manual washer for 12 x 75mm test tubes
- Produced from high-grade stainless steel and polypropylene,
 the washer is fully autoclavable as well as highly resistant to aggressive chemicals.
- Unique Two-in-One Design: Each tube consists of an inner tube that releases the distilled water washing solution into the wells for rinsing, while the outer tube vacuums away the liquid upon completion of the washing procedure
 - The vacuum pump is available for 115V and 230V connections
 - All components included in the kits can also be purchased individually
 - No programming required
 - No electrical wiring connection

Ordering Information

Cat. No.	Description
W-8	CAPPWash 8-channel
W-12	CAPPWash 12-channel
W-16	CAPPWash 16-channel
W-8KIT-115V	CAPPWash Kit including W-8, WP-115V, WB-1, WB-3, WB-4, and W-1000 (3 meters)
W-12KIT-230V	CAPPWash Kit including W-12, WP-230V, WB-1, WB-3, WB-4, and W-1000 (3 meters)

