

Single Channel Pipette Specifications

Color Code		Volume µL	Inaccuracy		Imprecision	
	Size		Absolute ± µL	Relative ± %	≤ SD, μL	≤ % CV
Violet	0.1-2µL	0.2 1 2	0.024 0.031 0.040	12.00 3.10 2.00	0.012 0.016 0.020	6.00 1.60 1.00
white	0.5-10µL	1 5 10	0.025 0.060 0.100	2.50 1.20 1.00	0.018 0.04 0.050	1.80 0.80 0.50
Grey	2-20µL	2 10 20	0.10 0.12 0.20	5.00 1.20 1.00	0.03 0.06 0.06	1.50 0.60 0.30
Red	5-50µL	5 25 50	0.10 0.20 0.50	2.00 0.80 1.00	0.06 0.09 0.15	1.20 0.35 0.30
Orange	10-100µL	10 50 100	0.25 0.45 0.70	2.50 0.90 0.70	0.10 0.15 0.20	1.00 0.30 0.20
Yellow	20-200µL	20 100 200	0.50 0.80 1.20	2.50 0.80 0.60	0.14 0.30 0.40	0.70 0.30 0.20
Blue	100- 1000µL	100 500 1000	2.00 4.00 8.00	2.00 0.80 0.80	0.60 0.75 2.00	0.60 0.15 0.20
Green	1-5mL	1000 2500 5000	12.00 20.00 30.00	1.20 0.80 0.60	3.00 5.00 7.50	0.30 0.20 0.15
Green	1-10mL	1000 5000 10000	30.00 40.00 60.00	3.00 0.80 0.60	6.00 10.00 15.00	0.60 0.20 0.15

Multichannel Pipette Specifications

Color Code	Size	Volume µL	Inaccuracy		Imprecision	
			Absolute ± µL	Relative ± %	≤ SD, µL	≤ % CV
Violet	0.2-2µL	0.2 1 2	0.048 0.062 0.080	24.00 6.20 4.00	0.024 0.032 0.040	12.00 3.20 2.00
White	0.5-10µL	1 5 10	0.050 0.116 0.200	5.00 2.32 2.00	0.036 0.064 0.100	3.60 1.28 1.00
Grey	2-20µL	2 10 20	0.10 0.23 0.40	5.00 2.30 2.00	0.072 0.128 0.20	3.60 1.28 1.00
Red	5-50µL	5 25 50	0.16 0.40 0.70	3.20 1.60 1.40	0.08 0.18 0.30	1.60 0.72 0.60
Orange	10-100µL	10 50 100	0.25 0.75 1.30	2.50 1.50 1.30	0.12 0.30 0.50	1.20 0.60 0.50
Yellow	20-200µL	20 100 200	0.36 1.44 2.40	1.80 1.44 1.20	0.16 0.44 0.80	0.80 0.44 0.40
Black	30-200µL	30 100 200	0.50 1.44 2.40	1.66 1.44 1.20	0.18 0.44 0.80	0.60 0.44 0.40
Black	125-300µL	125 150 300	1.76 2.10 4.20	1.40 1.40 1.40	0.55 0.54 0.90	0.44 0.36 0.30

Tested with distilled water at 19-22°C Test procedures in accordance to ISO 8655 All pipettes are tested with ExpellPlus" low retention tips, except 30-300 µL, 100-1000 µL, transferring of body heat. The transfer of heat affects 1-5mL and 1-10mL pipettes for which standard Expell tips are used.

Tips to improve the pipetting technique and results

Use the appropriate pipette tip size from a high quality manufacturer. CAPP pipettes are designed for optimal use with CAPP Expell pipette tips.

Work in a controlled environment. Environmental factors that affect pipetting volumes include: changes in temperature, changes in air pressure, vibrations and movement within the solution being pipetted.

Pre-rinsing tips helps the polypropylene tip, which is naturally hydrophobic, to adjust to the aqueous solution being pipetted. CAPP recommends three pre-rinse cycles of aspiration and dispense for each sample pipetted.

To eliminate unwanted liquid on the outside of the tip, it is recommended to touch-off the outside of the tip to the side of the vessel containing the sample. Unwanted droplets hanging from the bottom of the tip can be eliminated by touching the droplet to the surface of the aqueous solution.

Use consistent plunger pressure and speed for each sample.

Proper tip immersion depth is important for aspirating an accurate sample. Immersing the tip too deeply could cause extra liquid to be aspirated because of the additional force exerted by the aqueous sample. A shallow immersion depth could cause the tip to lose contact with the aqueous solution and aspirate a too small sample. The ideal tip immersion depth is between 2 to 5 mm during sample aspiration.

Immediately after aspirating the sample, pause for a second to allow the sample to fully equilibrate within the tip, before removing the tip from the solution.

Try to keep the pipette as vertical as possible during use. Do not incline the pipette more than 15° while aspirating or dispensing.

After dispensing the sample, touch-off the tip onto the wall of the receiving vessel container to ensure complete delivery of the sample.

Minimise handling of the pipette and tip to avoid the size of the air cavity in the tip and within the body of the pipette. It can affect the amount of sample aspirated and dispensed.

Pipettes Mechanical **JPP**

ecopipette

Recommended Volume Ranges

To maintain the highest level of accuracy and precision, it is important to use the appropriate instrument relative to the desired sample volume. Piston stroke length and the size of the air pocket in the tip are the two factors that make air displacement pipettes inaccurate when used at inappropriate volumes.

Piston Stroke Length:

Longer piston stroke lengths increase the accuracy and precision. A longer stroke length allows the error of inexact volumes to be dispersed over a larger volume range. Small errors are more impactful on a short stroke length than they are on a long stroke length. Piston stroke length is the reason that pipettes produce more accurate results on the high end of their volume range than they produce on the low end of their volume range.

Tip Air Pockets:

It is necessary to have an air pocket in a pipette tip. However, having too large of an air pocket will affect the accuracy of the results.

Fixed Volume Controller:

If you frequently pipette certain volumes, we recommended that you order additional fixed volume controllers to your specifications as this offers the safest and most convenient volume setting of all. Please list the desired volume(s) as well as the colour coding of the pipette when ordering.

General operation

Volume Adjustments (variable volume pipettes):

2) Choose the desired volume by turning the coloured

CAPP Expell tips are recommended for optimal use

Put the shaft into the tip and press firmly (Do not use

excessive force which will cause an over insertion of

the tip). Lightly tap the pipette tip back into the box

while loaded on the pipette shaft to ensure the tip

is properly seated and a good seal is formed. After dispensing the sample, eject the tip using the ejector

2) Place the tip in the liquid, release the plunger slowly

to aspirate and hold the tip in the liquid for 1 second.

3) Place the tip in the sample vessel. Push plunger to

4) Press the plunger to the second stop to blow out

5) Return the plunger back to the resting position.

3) Re-lock by pushing the lock lever upwards.

Loading and Ejecting Tips:

with the CAPP pipette.

Five Step Pipetting Cycle:

the first stop to dispense.

any remaining liquid.

Cleaning the Pipette:

Autoclaving:

1) Press the plunger down to the first stop.

downwards.

cap.

button.

Warranty or Service

The pipette is guaranteed for a period of three years from the date of delivery against defects in materials and workmanship of the product. The warranty covers the costs of defective materials and labour. The warranty does not cover common wear and tear or customer misuse. CAPP recommends servicing pipettes on an annual basis to keep them in optimal working condition.

All warranty work and repair must be done by a CAPP authorised service centre. To maintain the validity of the warranty, always have CAPP pipettes serviced by CAPP authorised service centres. The CAPP factory warranty is voided if repairs are performed by non-CAPP authorised service centre. A complete list of CAPP authorised service centres can be found at www.capp.dk.

Register Your CAPP Pipette

Your pipette has been engraved with a serial number (two letters and five digits) on the body of the pipette. Registering the pipette with CAPP will apply the warranty and create an ownership record for the instrument. Please refer to the serial number for all inquiries relating to the pipette.

Register Your Pipette!

Go to www.capp.dk and choose the pipette model you want to register. Click on 'Register your pipette'.

	Tro
 Endy sufficients investor 	dia
canada data canada da	F
Tara ta dattar	F
And an in the same is not the other	h
	F
 Nally participation through 	~
	Tang tig gjattan Tang tig gjattan Tang an tir datar gjattan filosoonid

Liquid is leaking from pipette tip	Pipette tip does not fit properly onto the pipette shaft	Use the appropriate size CAPP Exp	
	d is leaking from	Liquid being pipetted is hot or cold. Liquid being pipetted is very dense or viscous.	Shorten the amount of time the liquid Another possible solution is t use reverse mode pipetting
	Pipette shaft or sealing O-ring is worn	Replace shaft and/or sealing O-	
		Tip is not loaded properly to maintain a seal	Press on tip more firmly (avoid using force and over inserting the ti
Amount of sample delivered is not accurate	Pipette is not within calibration centre for specifications	Send the pipette to an authoris CAPP service centre cleaning and ma	
	Improper pipette technique	See the CAPP suggestions for imp pipetting technique and resul	
Uneven movement of piston or plunger	Piston has corrosion or contamination	Send the pipette to an authoris CAPP service centre for cleaning and m	
	Plunger rod has been damaged	Send the pipette to an authori CAPP service centre for cleaning and m	

1) Unlock the volume dial by pushing the lock lever

External cleaning is recommended using normal cleaning agents such as water with mild detergents, antiseptic agents or 60-70% ethanol. CAPP recommends that cleaning of internal components should be done by a trained professional in an authorised CAPP service centre.

Simply take off the volume controller and put the pipette in the autoclave. It must not exceed 121°C and a period of 20 minutes.

p pipettes. Our local star will contact pro to a free comple for trial.	
IN BOOMETR PRIMITER	
et piloverte kan kejt.	
B7-WoHtk	
Novembers	
BILLING COLORAD)

kpell tip

is in the tip. to

-ring

too much tip)

ised aintenance

proving ults

ised maintenance

ised maintenance







AHN Biotechnologie GmbH Uthleber Weg 14 | 99734 Nordhausen | Germany T: +49 (0) 3631 65242-0 | F: +49 (0) 3631 65242-90

www.capp.dk | info@cappahn.com