



CappOrigami™ Reagent Reservoirs

Fold, pour and pipette...!

- *V-shape ensures full usage of all reagents*
- *Flick-up design saves storage space*



The CappOrigami™ represents a innovative way of thinking when it comes to reagent reservoirs as subsequently described:

Perfect V-shape


The shape of the reservoir along with the stiffness of the material used ensures that the reservoir forms a sharp V-shape at the bottom. This makes your pipetting easier and also enables you to save on reagents as you only have to pour the amount needed into the reservoir.

Flat when stored


The CappOrigami™ reservoirs are packed as flat sheets. This way, they take up considerably less room in storage.

Easy to use


To use a CappOrigami™ reservoir, simply flick it up by pushing the sides together, which is very easy due to the prefolded lines. As soon as the CappOrigami™ reservoir reaches the proper shape you can pour liquid into the reservoir and start pipetting. The folds and the polypropylene material used assure that the reservoirs stay upright once you have folded them into the proper reservoir shape.

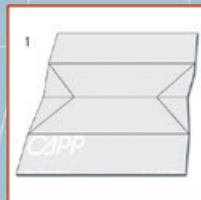

Take a flat CappOrigami™ reservoir sheet out of the pack.

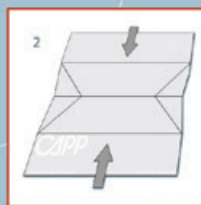

Push the two outer sides against the middle to flick the CappOrigami™ reservoir into shape.

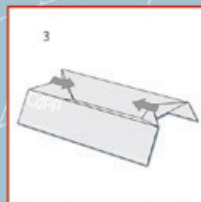


Push the end triangles in to make the CappOrigami™ reservoir ready for use.

We take pipetting to perfection.

Ordering information:

Code	Description	Unit	Packaging
CA40505	30 ml reservoirs for 8- and 16-channel pipettes	Pack/50	Non-Sterile
CA40506	30 ml reservoirs for 8- and 16-channel pipettes	Pack/5x10	Sterile
CA40510	40 ml reservoirs for 12-channel pipettes	Pack/50	Non-Sterile
CA40511	40 ml reservoirs for 12-channel pipettes	Pack/5x10	Sterile

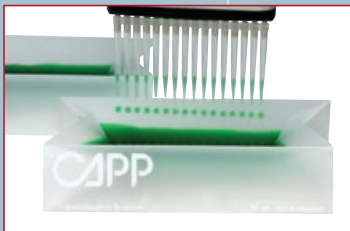

CappAero™ Multi

Standard	Type	Volume	Colour
C10-8	8-channel	0.5-10 µl	○ White
C20-8	8-channel	2-20 µl	● Grey
C50-8	8-channel	5-50 µl	● Red
C100-8	8-channel	10-100 µl	● Orange
C200-8	8-channel	20-200 µl	● Yellow
C300-8	8-channel	30-300 µl	● Black
C10-12	12-channel	0.5-10 µl	○ White
C20-12	12-channel	2-20 µl	● Grey
C50-12	12-channel	5-50 µl	● Red
C100-12	12-channel	10-100 µl	● Orange
C200-12	12-channel	20-200 µl	● Yellow
C300-12	12-channel	30-300 µl	● Black
C02-16	16-channel	0.2-2 µl	● Violet
C10-16	16-channel	0.5-10 µl	○ White
C50-16	16-channel	5-50 µl	● Red



CAPP®

Capp A/S

Landbrugsvej 10 • DK 5260 Odense S
Denmark
tel. +45 6613 6140 • fax. +45 6613 2770
www.capp.dk • info@capp.dk

Distributor